



Uniblitz[®] FS6

6mm Uni-Stable Laser Shutter

Overview

The Uniblitz FS series shutters are designed and optimized to operate directly from +5VDC and do not require a separate driver. Removing the +5VDC (0VDC) closes these shutters. A simple control circuit can be used to operate these shutter devices from a TTL trigger pulse. This control can be also accomplished with our new **VLM1 TTL control interface module**, which is now available.

This low-cost innovation provides the reliability of Uniblitz shutters (typical lifetime >1 million cycles) at a single operating voltage.

Key Features

- 6mm aperture
- Default closed operation
- RoHS Compliant
- Transfer time on opening:3.3 milliseconds
- Transfer time on closing:6.0 milliseconds

Specifications

Electrical Specifications	
Coil resistance	12 OHMS
Voltage to Open	+5 VDC at 425 mA
Hold Voltage	+5 VDC

Mechanical Specifications					
Weight Unhoused	25.7 g				
Weight Half-Housed	59.5 g				
Weight Housed	94.5 g				
Operating Temp.	-40 - +65 C				
Max. Opening Bounce	15%				
Max. Closing Bounce	5%				
Max. Freq. of Operation ¹	50 Hz				
Number of Shutter Blades	1				

¹ (Continuous/Burst) Continuous frequency rating specified at shutter's minimum exposure time (MET).



Need Support? Please visit our website or email us at **info@uniblitz.com**. Tel: +1 585-385-5930 | Fax: 585-385-6004 | 803 Linden Ave. Rochester, NY 14625 Updated 8/23 | Document Version 6.4 | ©2023 Vincent Associates

Shutter Timing Data



MET¹ and TET² for FS6

- Minimum Exposure Time (MET): 10.0 msec. ³
- Typical Exposure Time (TET):
 20.0 msec. ³

FS6 4 (w/ VLM1 (+5 VDC) and "T" blades) 5		Min. Exp. Time	Min. Exp. Time (msec.)		Typ. Exp. Time (msec.)	
Graph	Description	Mean	Max	Mean	Мах	
0 - A	Delay time on opening after current is applied	4.5	5.5	4.5	5.5	
A - C	Transfer time on opening	3.5	4.5	3.5	4.5	
0 - C	Total opening time	8.0	10.0	8.0	10.0	
B - F	Minimum equivalent exposure time	7.7	9.5	19.3	21.0	
С - Е	Minimum dwell time (based on exposure pulse)	4.0	4.5	15.5	16.0	
D - E	Delay time on closing after current is removed	2.0	4.5	3.5	6.0	
E - G	Transfer time on closing	3.8	5.5	4.0	5.5	
A - G	Total window time	11.3	14.5	23.0	26.0	

¹ The minimum exposure pulse required to achieve a full opening and closing.

² The exposure pulse required to ensure that the A - G time will change linearly with the exposure pulse (E.g. 1 msec changes in exposure pulse beyond 20 msec will provide 1 msec changes in A-G time).

³ Exposure pulses provided between 10 msec and 20 msec will open and close the shutter fully, however timing changes between input and output may not be linear. This is due to shutter design, construction, and adjustment.

⁴ Individual timing segments may vary.

⁵ Under no circumstances should any type of lubricant be applied to the shutter blade area. Lubricating the shutter blades will likely slow the shutter down and may eventually render it inoperable.

Uniblitz® FS6

Product Options

FS6 **23456**-7

Ex: FS6S2Z0L-EC



² Various mounting methods available depending on housing option.

³ Other blade coating options may be available by special order.

⁴ Input side only; Teflon® coating is on opposite side to protect shutter blade surface. Light source must be input to the reflective side only.

⁵ Vacuum compatibility up to 10E-10 Torr.

Electronic Sync.

The synchronization system for FS shutter devices incorporates a small magnet mounted to the driving mechanism and a Hall effect sensor. When the device achieves approximately 80% of full open, the magnet causes the Hall effect sensor to change state, producing a signal to indicate that the shutter has switched to the active state. Shown to the left is the FS series shutter schematic which incorporates the electronic synchronization system. There is no connection to the designated synchronization pins when an electronic sync. is not selected.



Typical Control Circuit

This simple control circuit to the right can be used to operate the shutter device from a TTL trigger pulse.¹ This control can also be accomplished with the **VLM1 interface module.**

¹ Note that the shutter will not operate directly from a TTL signal.



Half-Housing - FS6S2T0L



Un-Housed - FS6S1T0L









Uniblitz® FS6 Technical Drawings

Full-Housing - FS6S3T0L



- OPTIONAL REFLECTIVE SURFACE OPPOSITE
- ACTUATOR COIL SIDE.