



Uniblitz[®] VS35

35mm Uni-Stable Optical Shutter

Overview

The Uniblitz VS35 has been designed to provide accurate, repeatable exposures for a wide variety of applications. The VS series shutters have been successfully field tested in microscopy, video imaging, PMT protection, and photographic applications worldwide. The VS35 is available in a housed or unhoused configuration for OEM applications (or where spatial limitations are a consideration). Uni-stable shutter devices, like the VS35, require power to hold the blades in the open state (though the device can be custom-configured for "normally open" operation).

Key Features

- 35mm aperture
- Small, slim form-factor
- Configured for the <u>VCM-D1</u>
 <u>Shutter Driver</u>
- RoHS Compliant
- Transfer time on opening:
 13.0 milliseconds
- Transfer time on closing:
 12.0 milliseconds

Specifications

Electrical Specifications	
Coil resistance	12 OHMS
Voltage to Open	+70 VDC
Hold Voltage (Nominal) ¹	+7 VDC / +5 VDC ²

Mechanical Specifications	
Weight Unhoused	70.0 g
Weight Housed	410.0 g
Operating Temp.	0 - 80 °C
Max. Opening Bounce	15%
Max. Closing Bounce	5%
Max. Freq. of Operation ³	5 Hz / 20 Hz
Number of Shutter Blades	2

¹*Voltage level required across actuator coil when being held in the open position.*

² Dual hold voltage system included in the VCM-D1 Driver.

³ (Continuous/Burst) Continuous frequency rating specified at shutter's minimum exposure pulse. Burst frequency rating specified for four (4) seconds maximum with one (1) minute minimum between bursts.



Need Support? Please visit our website or email us at **info@uniblitz.com**. Tel: +1 585-385-5930 | Fax: 585-385-6004 | 803 Linden Ave. Rochester, NY 14625 Updated 5/24 | Document Version 6.0 | ©2024 Vincent Associates

Uniblitz[®] VS35

Shutter Timing Data



¹ Under no circumstances should any type of lubricant be applied to the shutter blade area. Lubricating the shutter blades will likely slow the shutter down and may eventually render it inoperable.

VS35 (w/ VCM-D1 and "T" blades) ¹ Time	e (msec.)
0 - A	Delay time on opening after current applied	d 5.0
A - C	Transfer time on opening	13.0
0 - C	Total opening time	18.0
C - E	Min. dwell time with min. input pulse	5.0
B - F	Min. eqivalent exp. time	20.0
E - G	Transfer time on closing	12.0
A - G	Total window time	30.0
MET	Min. exposure time	20.0
TEP	Typical exposure pulse	>23.0

Product Options

VS35 23456-7-8

Ex: VS35S2ZM0-EC-21



² Other blade coating options may be available by special order.

³ Input side only; Teflon[®] coating is on opposite side to protect shutter blade surface. Light source must be input to the reflective side only.

⁴ Mounts are only compatible if #2 housing is optioned as well.

Uniblitz[®] VS35

Electronic Sync.

The electronic synchronization system provides a feedback signal (through the driver utilized) after the shutter transfers to the open state. The system incorporates an infrared emitting diode, an infrared sensitive detecting transistor, and an interrupting vane. The vane is attached to the shutter so as to block the light path between the emitter and detector in the closed position. When the shutter transfers to the 80% open position, the vane is removed from the infrared light path, allowing the emitter to switch the detector to the active state. For the VS35, this system uses a similar activation flag attached to the mechanism, which triggers a reflective emitter/ detector device. No connection to the designated synchronization pins when no electronic sync. is selected.



Uniblitz® VS35 Technical Drawings

Un-Housed



Housed



- (3) #2-56 MOUNTING HOLES ON A Ø 1.578 BOLT CIRCLE, 120° APART
- (3) #4-40 MOUNTING HOLES ON A Ø2.425 BOLT CIRCLE 120° APART (UNDER NAME PLATE)
- ALL DIMENSIONS [MM] INCH





Uniblitz[®] VS35 Technical Drawings

Connector Layout





VIEW B-B

NOTES:

- 1. DRAWING SHOWING CONNECTOR AND 1/4-20 MOUNTING LOCATION.
- 2. FINISH BLACK ANODIZE.
- 3. TERMINATES TO 710P OR 710C CABLE ASSEMBLY.
- 4. ALL DIMENSIONS MAXIMUM [MM]

